Consort

MANUAL

HANDLEIDING

MODE D'EMPLOI

ANLEITUNG

C6010
C6020
C6030

Table of contents

Introduction	1
Manufacturer	1
Warranty	1
Servicing.....	1
Keyboard	1
Language	2
Outputs	2
Contrast	2
Date and time	2
Password	2
User standard tables	2
mV measurement	3
Temperature measurement	3
pH measurement.....	3
Important	3
Conductivity measurement.....	4
Dissolved oxygen measurement	5
Principle.....	5
Interferences	5
Maintenance of pH electrodes.....	6
Maintenance of metal electrodes	6
Maintenance of conductivity electrodes	6
Maintenance of oxygen electrodes.....	6
Ion measurement	7
GLP.....	7
Calibration reminder	7
Data-logging	8
USB.....	8

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Introduction

This instrument is manufactured with the latest technology and needs no particular maintenance. **Consort** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

The normal operating temperature should be between 0° and 40°C. Never use the instrument in a room with high humidity (>95 %) or at very low temperatures (condensation water!).

Manufacturer

Consort bvba

Parklaan 36

B2300 Turnhout

Belgium

Tel (+32)(14)41 12 79

Fax (+32)(14)42 91 79

E-mail: info@Consort.be

Warranty

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **Consort** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **Consort** before returning any product for warranty repair on a freight prepaid basis!

Consort is not liable for consequential damages arising out of the use or handling of its products.

Servicing

In the event of this instrument being returned for servicing, the owner is requested to remove the power supply lead and **NOT** to send the following items unless they are suspect:

Manual

Cables

Accessories

If serious malfunctioning occurs, stop using the unit immediately and consult your local **Consort** dealer.

Batteries

The batteries can be replaced by opening the bottom compartment of the cabinet. Only use NiMH batteries!

AC adaptor

Connect the AC adaptor (9 V, 300 mA for 230 V~ or 120 V~) to the DC socket. **Avoid to use an AC adaptor without NiMH batteries being inserted in the instrument!** Do not hold the adaptor by wet hand!

Keyboard

MODE = Selects the settings or escapes from error traps, calibration procedures, etc..by returning to the original mode.
↑↓ = Buttons for entering a value or for selecting a function.
OK = Proceeds a function.
CAL = Starts a calibration.
STORE = Stores the displayed value or send it to a computer or printer.
HOLD = Holds display when measuring or shows a built-in manual during a setup.

The measuring electrode should be connected to the coaxial INPUT connector. If separate electrodes are used, connect the reference electrode to the REF. terminal. Automatic temperature compensation and temperature measurements are possible by plugging a Pt1000 temperature probe into the °C terminals. You can also use a combination pH electrode with built-in Pt1000. Its banana plugs should be inserted in the °C terminals. Without Pt1000, the manual temperature compensation is automatically switched on.

An insulated standard USB port is provided for interfacing the instrument with a computer.

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [SYSTEM] and press **OK**.
4. Select [LANGUAGE] and press **OK**.
5. Select the desired language and press **OK**.

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [SYSTEM] and press **OK**.
4. Select [CONTRAST] and press **OK**.
5. Adjust the contrast and press **OK**.

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [SYSTEM] and press **OK**.
4. Select [DATE/TIME] and press **OK**.
5. Select [MODIFY] and press **OK**.
6. Adjust date and time and press **OK**.

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [SYSTEM] and press **OK**.
4. Select [PASSWORD] and press **OK**.
5. A private code can be programmed to avoid undesired access to the instrument. Select [YES] and press **OK** to enter your secret sequence of 5 keys.

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [SYSTEM] and press **OK**.
4. Select [POWER MANAGEMENT] and press **OK**.
5. Select [POWER OFF] (batteries or DC power) and press **OK**.
6. Adjust the time to shut down and press **OK**.
7. Select [BACKLIGHT] (only in case of DC power) and press **OK**.
8. Select [YES] or [NO] and press **OK**.

Important

- A blinking decimal point warns you for unstable measurements. Wait to read the display!
- Stirring the solution during the measurements promotes the homogeneity and is obligatory!
- The instrument will refuse automatic calibration when the electrode is unstable. Insufficient stirring or a worn electrode may be the cause.

pH measurement

1. Select the desired mode (pH) by pressing **MODE**. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
2. The instrument permits to choose between the buffers in memory (1.68, 2.00, 4.00, 4.01, 6.87, 7.00, 9.18, 9.21, 10.01, 12.00, 12.45 + max. 5 user tables). Select the proper values and press **OK**. The unused buffers should be switched off.
3. Rinse the electrodes with distilled water and immerse them in the first buffer solution. Select [**CALIBRATE**], press **OK** and follow the instructions on the screen until the calibration is finished.
4. After rinsing the electrodes with distilled water, immerse them in the samples and read the display.
5. Rinse the electrodes always with distilled water after use and store them in a 3...4 M KCl solution.

Using pH electrodes with an abnormal zero point:

1. Press **CAL**.
 2. Select [**Eo = x mV**] and press **OK**.
 3. Enter the zero point of the electrode and press **OK**.
- *Hold **MODE** pressed and press ↓ to change the resolution from 0.1 to 0.001 pH.*

mV measurement

1. Select the desired mode (mV) by pressing **MODE**. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
 2. Immerse the electrodes in a standard solution of known potential. Select [**CALIBRATE**] and press **OK**. Adjust to the proper value and press **OK**.
- *Select [**RESET**] and press **OK** to reset the calibration.*
 - *Hold **MODE** pressed and press ↓ to change the resolution from 1 to 0.1 mV.*

Temperature measurement

1. Select the desired mode (°C) by pressing **MODE**. Should you want to recalibrate, press **CAL**.
 2. Immerse the Pt1000 in a solution of known temperature. Select [**CALIBRATE**] and press **OK**. Adjust to the proper value and press **OK**.
- *Select [**RESET**] and press **OK** to reset the calibration.*
 - *Without Pt1000, press **CAL**, adjust the manual temperature compensation and proceed by pressing **OK**.*

1. Select the desired mode (S/cm) by pressing **MODE**. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
 2. The instrument permits to choose between the standards in memory (1413 $\mu\text{S/cm}$, 12.88 mS/cm, 111.8 mS/cm + max. 3 user tables). Select the proper values and press **OK**. The unused standards should be switched off.
 3. After rinsing the electrode several times with the first standard solution, immerse it in that standard. The solution temperature is not so critical but should lie between 0°C and 30°C. When no Pt1000 is used, do not forget to compensate manually first! Select [CALIBRATE], press **OK** and follow the instructions on the screen until the calibration is finished.
 4. Rinse the electrode several times with the sample, immerse it in that solution and read the display.
 5. Rinse the electrode always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).
- *When starting a titration, hold **MODE** pressed and press \downarrow to lock the actual range and avoid cross-over errors due to differences in the measuring frequency. Hold **MODE** pressed and press \downarrow to return again to the normal measuring mode.*

Temperature compensation:

1. Press **CAL**.
2. Select [REF.] and press **OK**.
3. Select the temperature to which all future measurements will be referred to (25°C or 20°C) and press **OK**.

Capacitive compensation:

1. Capacitive compensation increases accuracy in the very low conductivity ranges (<10 $\mu\text{S/cm}$). Select whether or not this compensation should be applied. Verify if the attached electrode is completely dry and press **CAL**.
2. Select [CAP.COMP] and press **OK**.
3. Select [YES] or [NO] and press **OK**. Follow the instructions on the screen.

Resistivity measurement:

1. Select the desired mode ($\Omega\cdot\text{cm}$) by pressing **MODE**. The display will immediately show the measured value.

Salinity measurement:

1. Select the desired mode (SAL) by pressing **MODE**. The display will immediately show the measured value according to the Practical Salinity Scale (PSS-78) which is based upon standard seawater at 15°C.

TDS measurement:

1. Select the desired mode (mg/l)(TDS) by pressing **MODE**. The display will immediately show the measured value.
- *To browse between conductivity, resistivity, salinity or TDS, hold **MODE** pressed and press \uparrow sequentially.*

Dissolved oxygen measurement

(C6020, C6030)

1. Select the desired mode (ppm) or (%O₂) by pressing **MODE**. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
 2. The display shows the salinity correction. Select the proper value and press **OK**. Leave salinity correction to zero unless you are going to measure in heavily salted solutions such as e.g. sea-water (35 g/l). Select [CALIBRATE], press **OK** and follow the instructions on the screen until the calibration is finished.
 3. The electrode exposed to the air reaches an equilibrium corresponding to the partial pressure of oxygen and thus to saturation in water at the given temperature. The instrument shows the measured saturation, current, temperature, and will calibrate automatically when readings are stable.
 4. After rinsing the electrodes with distilled water, immerse them in the samples and read the display. Stirring the solution during the measurements promotes the homogeneity and is **obligatory!** The advection rate must be at least 10 cm/s.
 5. Rinse the electrode always after use and store it in distilled water.
- *A blinking decimal point warns you for unstable measurements. Wait to read the display!*
 - *Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.*
 - *Hold **MODE** pressed and press ↓ to change the resolution from 0.1 to 0.01 ppm or from 1 to 0.1 %.*

Air pressure measurement:

1. Select the desired mode (hPa) by pressing **MODE**. Should you want to recalibrate, press **CAL**.
 2. Select [CALIBRATE], press **CAL**, adjust to the real air pressure and press **OK**.
- *Select [RESET] and press **OK** to reset the calibration.*

Principle

(C6020, C6030)

The oxygen meter and its electrode function according to the Clark principle with silver as cathode and lead as anode in an electrolyte cell. Oxygen gas present in the electrolyte is reduced to OH ions at the cathode. The resulting current is diffusion limited and therefore proportional to the oxygen concentration in the sample solution. This current is amplified, corrected, and displayed in mg/l, ppm or % dissolved oxygen.

Interferences

(C6020, C6030)

All substances which can diffuse through the membrane and for which 800 mV potential suffices for polarographic reduction, will be reduced in the electrode. This will give a corresponding current contribution, if they are present. Interference can be caused by ions entering the electrode through porous or mechanically damaged membranes and by diffusion of other reactive gases apart from oxygen, e.g. CO₂, Cl₂, SO₂, and H₂S. These substances react in undesired manner with the electrode. Acidic or basic gases change the pH value of the electrolyte solution and thus disturb the reading, particularly when measuring small oxygen concentrations. High salt concentrations in the sample solution can falsify readings too.

A pH electrode is active and stable only after wetting! For this purpose it must be immersed for **at least ten hours** in a 3...4 M KCl solution. During short interruptions (e.g. storage) the electrode should be immersed in a 3...4 M KCl solution. In doing this it is always kept ready for use. When the interruption is longer than a month, refill the closing cap with 3...4 M KCl and plug it on the electrode tip in order to protect the glass bulb. Before use, ensure that the reference part of the electrode is topped up with a 3...4 M KCl solution.

Avoid a low pressure inside the electrode! Therefore always remove the closure from the refilling aperture during the measurements as well as during the calibration. This allows the saltbridge solution to flow through the ceramic liquid junction and prevents contamination of the electrolyte. For the same reason, the inside level should always be higher than the outside level of the measuring solution. Close the refilling aperture again when storing the electrode.

A polluted electrode may be cleaned with a soft detergent or 0.1 M HCl. Greasy substances may be removed with acetone or alcohol (**never do this with plastic electrodes!**).

If the electrode is polluted by proteinaceous materials (such as blood), it should stand in a cleaning solution overnight and then be cleaned with distilled water before use. The pH electrode wears away by being used. If the electrode tends to respond slower and calibration becomes difficult, even after cleaning, it should be replaced by a new one.

Maintenance of metal electrodes

Metal electrodes (Pt, Ag, Au): Metal electrodes are always ready for use. During short interruptions they are immersed in distilled water. **They should be cleaned regularly:**

- Silver electrodes are immersed in a concentrated ammonia solution during one hour.
- Platinum or gold electrodes are immersed in concentrated nitric acid during one hour.

Maintenance of conductivity electrodes

A conductivity cell is active and stable only after wetting! For this purpose it must be immersed for **at least one hour** in distilled water. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

A polluted cell may be cleaned with a soft detergent or diluted nitric acid. Greasy substances may be removed with acetone or alcohol (never do this with plastic electrodes!).

Maintenance of oxygen electrodes

A dissolved oxygen electrode is active and stable only when polarised! The electrode body has almost unlimited life and requires no maintenance. After prolonged use of the electrode, it may become deactivated. An indication is that the electrode no longer responds correctly to calibration. In this case:

1. Carefully screw off the measuring head.
 2. Cautiously remove the precipitated deposits with the aid of filter paper. **Do not use grinding paper or a glass fibre!**
 3. Rinse several times with distilled water and shake carefully off the water drops.
 4. Take the measuring head, fill it with new electrolyte solution and very slowly screw it onto the electrode while holding it vertically.
 5. **Let the electrode rest for a few hours!** The electrode is now ready for use again.
- *Replace the membrane only when damaged, not when calibration is no longer possible!*

Ion measurement

(C6030)

1. Select the desired mode by pressing **MODE**. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
 2. The instrument permits to choose between the possible standards (1, 10, 100 ng/l, 1, 10, 100 µg/l, 1, 10, 100 mg/l, 1, 10, 100 g/l). Select the proper values and press **OK**. The unused standards should be switched off.
 3. Select [**CALIBRATE**] and press **OK**.
 4. After rinsing the electrodes with distilled water, immerse them in the first standard solution. Select the first calibration value and press **OK** when readings are stable. Go on in the same way with the next standard solutions.
 5. Decide whether a blank correction should be carried out or not. Follow the instructions on the screen. After rinsing the electrodes with distilled water, immerse them in a blank solution. When readings are stable press **OK**.
 6. Rinse the electrodes with distilled water, immerse them in the samples, and read the concentration on the display.
- *Maintenance: we refer to the manual supplied with the ion specific electrodes.*

GLP

1. Select the desired range by pressing **MODE** and then press **CAL**.
2. Select [**GLP**] and press **OK**.
3. Select [**SHOW REPORT**] and press **OK**. Browse with **↑↓** to show a complete calibration report.
Select [**SEND REPORT**] and press **OK** to send the report to a computer.

Calibration reminder

1. Select the desired range by pressing **MODE** and then press **CAL**.
2. Select [**GLP**] and press **OK**.
3. Select [**INTERVAL**] and press **OK**.
4. Select the desired time interval between each automatic warning for a new calibration of the electrodes and press **OK**.

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [DATA] and press **OK**.

Start the data-logging:

1. Select [ACTIVATE] and press **OK**.
2. Decide whether data-logging should be carried out or not and press **OK**.
3. Select [CONTINUOUS] and press **OK**.
4. Decide whether the oldest data should be overwritten when the memory limit is reached or not and press **OK**.
5. Select [INTERVAL] and press **OK**.
6. Select the desired time interval between the data-logging and press **OK**.
7. Return to the measurements by pressing **MODE**.
8. The data-logging starts according to the previous settings while the display shows the logging-number e.g. [LOG.00027].

Process the stored values:

1. Select [PROCESS] and press **OK**.
2. Select [TABLE] or [SEND] to display or send the stored data and press **OK** to continue. Follow the instructions on the screen.

Erase the stored values:

1. Select [PROCESS] and press **OK**.
2. Select [ERASE] and press **OK**. Follow the instructions on the screen.

USB

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [USB] and press **OK**.
4. Select [INTERVAL] and press **OK**.
5. Select the desired interval between the transmitted data and press **OK**. Pre-set to zero if no automatic transmitting is required.
6. Select [BAUD] and press **OK**.
7. Select the desired baudrate and press **OK**.

- *You can download a suitable USB driver from www.consort.be.*

User standard tables

1. Press **MODE**.
2. Select [MENU] and press **OK**.
3. Select [USER TABLES] and press **OK**.
4. Select [pH buffers] or [S/cm standards] and press **OK**.
5. Follow the instructions on the screen to enter the different values. The value for 25°C is obligatory!

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Specifications		C6010	C6020	C6030
pH	Range	-2...+16 pH	-2...+16 pH	-2...+16 pH
	Resolution	0.001 pH	0.001 pH	0.001 pH
	Accuracy	0.1% ± 1 digit	0.1% ± 1 digit	0.1% ± 1 digit
	Calibration	1...5 points	1...5 points	1...5 points
	Buffers	9 pre-programmed 5 user specified	9 pre-programmed 5 user specified	9 pre-programmed 5 user specified
	Temperature compensation	-5...+105 °C	-5...+105 °C	-5...+105 °C
	ISO-pH	6...8 pH	6...8 pH	6...8 pH
	Slope	80...120%	80...120%	80...120%
	Zero point (Eo)	±999 mV	±999 mV	±999 mV
mV	Range	±2000 mV	±2000 mV	±2000 mV
	Resolution	0.1 mV	0.1 mV	0.1 mV
	Accuracy	0.1% ± 1 digit	0.1% ± 1 digit	0.1% ± 1 digit
	Calibration	1 point	1 point	1 point
ION	Range			0.01 ng/L...100 g/l
	Resolution			3 digits
	Accuracy			0.5% ± 1 digit
	Calibration			2 points + blank
CONDUCTIVITY	Range (cc dependent)	0...2000 mS/cm	0...2000 mS/cm	0...2000 mS/cm
	Resolution (cc dependent)	0.001 µS/cm	0.001 µS/cm	0.001 µS/cm
	Accuracy	0.5% f.s. of range	0.5% f.s. of range	0.5% f.s. of range
	Calibration	1...3 points	1...3 points	1...3 points
	Standards	0.01/0.1/1M KCl 3 user specified	0.01/0.1/1M KCl 3 user specified	0.01/0.1/1M KCl 3 user specified
	Cell constant (cc)	0.1/1/10 cm ⁻¹ ±30%	0.1/1/10 cm ⁻¹ ±30%	0.1/1/10 cm ⁻¹ ±30%
	Temperature compensation	-5...+105 °C	-5...+105 °C	-5...+105 °C
	Reference temperature	20° or 25 °C	20° or 25 °C	20° or 25 °C
	Temperature coefficient	natural waters (EN27888)	natural waters (EN27888)	natural waters (EN27888)
	Range lock	✓	✓	✓
	Capacitive compensation	✓	✓	✓
	RESISTIVITY	Range	0...200 MΩ.cm	0...200 MΩ.cm
Resolution		1 Ω.cm	1 Ω.cm	1 Ω.cm
SALINITY	Range	0...70 ppt	0...70 ppt	0...70 ppt
	Resolution	0.1 ppt	0.1 ppt	0.1 ppt
TDS	Range	0...100 g/l	0...100 g/l	0...100 g/l
	Resolution	0.01 mg/l	0.01 mg/l	0.01 mg/l
DISSOLVED OXYGEN	Range		0...60 mg/L (0...600%)	0...60 mg/L (0...600%)
	Resolution		0.01 mg/L (0.1%)	0.01 mg/L (0.1%)
	Accuracy		1% ± 1 digit	1% ± 1 digit
	Calibration		1 point	1 point
	Temperature compensation		0...50 °C	0...50 °C
	Salinity compensation		0...40 ppt	0...40 ppt
TEMPERATURE	Range	-5...+105 °C	-5...+105 °C	-5...+105 °C
	Resolution	0.1 °C	0.1 °C	0.1 °C
	Accuracy	0.3 °C	0.3 °C	0.3 °C
	Calibration	1 point	1 point	1 point
AIR PRESSURE	Range		600...1300 hPa	600...1300 hPa
	Calibration		1 point	1 point
INPUTS	Measurement	BNC, 10 ¹² Ω	BNC, 10 ¹² Ω	BNC, 10 ¹² Ω
	Temperature	2 banana, for Pt1000	2 banana, for Pt1000	2 banana, for Pt1000
CALIBRATION	Reminder	0...999 h	0...999 h	0...999 h
	GLP	✓	✓	✓
DISPLAY	LCD	128x64 pixels	128x64 pixels	128x64 pixels
	White backlight	✓	✓	✓
	Hold function	✓	✓	✓
	Selectable resolution	✓	✓	✓
	Real time clock	✓	✓	✓
COMMUNICATION	Interface with computer	USB, insulated	USB, insulated	USB, insulated
DATA-LOGGING	Data sets	12000	12000	12000
	Modes	all (9)	all (9)	all (9)
	Manual or timed	✓	✓	✓
	Interval	1...9999 s	1...9999 s	1...9999 s
SECURITY	Identification number	✓	✓	✓
	Password protection	✓	✓	✓
AMBIENT CONDITIONS	Temperature	0...40 °C	0...40 °C	0...40 °C
	Humidity	0...95%, non condensing	0...95%, non condensing	0...95%, non condensing
POWER SUPPLY	Mains	210...250 VAC, 50/60 Hz	210...250 VAC, 50/60 Hz	210...250 VAC, 50/60 Hz
	Low voltage	9...15 VDC	9...15 VDC	9...15 VDC
	Batteries	4x1.2 V, AA, NiMH	4x1.2 V, AA, NiMH	4x1.2 V, AA, NiMH
DIMENSIONS	WxDxH	12x25x5 cm	12x25x5 cm	12x25x5 cm
WEIGHT	Meter	600 g	600 g	600 g

Temp.	1.68	4.01	6.87	9.18	12.45	<i>NIST buffers</i>	
0° C	1.666	4.003	6.984	9.464	13.423	(pH)	
5° C	1.668	3.999	6.951	9.395	13.207		
10° C	1.670	3.998	6.923	9.332	13.003		
15° C	1.672	3.999	6.900	9.276	12.810		
20° C	1.675	4.002	6.881	9.225	12.627		
25° C	1.679	4.008	6.865	9.180	12.454		
30° C	1.683	4.015	6.853	9.139	12.289		
40° C	1.694	4.035	6.838	9.068	11.984		
50° C	1.707	4.060	6.833	9.011	11.705		
60° C	1.723	4.091	6.836	8.962	11.449		
70° C	1.743	4.126	6.845	8.921			
80° C	1.766	4.164	6.859	8.885			

Temp.	2.00	4.00	7.00	9.21	10.01	12.00	<i>Standard buffers</i>	
5° C	1.85	3.99	7.08	9.45	10.24	12.56	(pH)	
10° C	1.89	3.99	7.06	9.38	10.18	12.35		
20° C	1.96	3.99	7.02	9.26	10.06	12.14		
25° C	2.00	4.00	7.00	9.21	10.01	12.00		
30° C	2.03	4.01	6.99	9.16	9.87	11.89		
40° C	2.08	4.03	6.98	9.06	9.89	11.71		
50° C	2.12	4.06	6.97	8.99	9.83	11.52		
60° C	2.14	4.09	6.98	8.93	9.79	11.33		

Temp.	0.01 M	0.1 M	1 M	<i>Conductivity standards</i>			
0° C	0.776	7.15	65.4	(mS/cm)			
5° C	0.896	8.22	74.1				
10° C	1.020	9.33	83.2				
15° C	1.147	10.48	92.5				
20° C	1.278	11.67	102.1				
25° C	1.413	12.88	111.8				
30° C	1.552	14.12	121.7				

Temp.	124	358	<i>Redox standards</i>				
0° C	156	388	(mV)				
5° C	148	381					
10° C	141	376					
15° C	135	370					
20° C	130	364					
25° C	124	358					
30° C	122	352					
35° C	116	346					
40° C	99	340					

DECLARATION OF CONFORMITY

We declare under our sole responsibility that the product

**Multi-channel analysers
content of the type numbers**

C6010, C6020, C6030

to which this declaration relates is in conformity
with the following standards

EN61010

LOW VOLTAGE DIRECTIVE 73/23/EEG

EN50081-1

EN50082-1

EN60555-2

EMC DIRECTIVE 89/336/EEG

Consort bvba

Parklaan 36, B-2300 Turnhout, Belgium

Tel: (+32) (0)14 41 12 79

Fax: (+32) (0)14 42 91 79

Sales: sales@consort.be

The electro-magnetic susceptibility has been chosen at a level that gains proper operation in residential areas, on business and light industrial premises and on small-scale enterprises, inside as well as outside of buildings. All places of operation are characterised by their connection to the public low voltage power supply system.

WARRANTY CERTIFICATE

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. Accessories and breakable items such as electrodes are not warranted unless proven to be defective before shipment.

The original purchase order numbers, Consort invoice numbers and serial numbers of the products must be provided.

CONSORT will repair all defective equipment returned to it during the warranty period without charge (CIF Turnhout prepaid by sender), provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures described in the instruction manual and when not having been subject to accident, alteration, misuse or abuse.

If the products have been used with or have come into contact with fluids, an MSDS (material safety data sheet) must be supplied prior to issuing a return authorisation.

A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

Consort

www.consort.be

Parklaan 36 • B-2300 Turnhout • Belgium

Tel: (+32) (0)14 41 12 79

Fax: (+32) (0)14 42 91 79

Sales: sales@consort.be

Support: support@consort.be

Information: info@consort.be